

IdeZ Protease (IgG-specific)



P0770S



4,000 units Lot: 0031512 Exp: 12/17
80,000 U/ml Store at -20°C

Description: IdeZ Protease (IgG-specific) is a recombinant antibody specific protease cloned from *Streptococcus equi* subspecies *zooepidemicus* that recognizes all human, sheep, monkey, and rabbit IgG subclasses, specifically cleaving at a single recognition site below the hinge region, yielding a homogenous pool of F(ab')_2 and Fc fragments. IdeZ Protease more effectively cleaves murine IgG2a than IdeS.

Cleavage Sites:

human IgG1, IgG3, IgG4: CPAPPELLG ∇ GPSVF
human IgG2: CPAPPVA ∇ GPSVF
murine IgG2a: CPAPNLLG ∇ GPSVF
murine IgG3: CPPGNILG ∇ GPSVF


Source: Cloned from *Streptococcus equi* subspecies *zooepidemicus* and expressed in *E. coli*.

Supplied in: 20 mM Tris-HCl, 50 mM NaCl,
1 mM EDTA, pH 7.5 @ 25°C

Reagents Supplied with Enzyme:
10X GlycoBuffer 2

Unit Definition: One unit is defined as the amount of enzyme required to cleave > 95% of 1 μg of human IgG, in 15 minutes at 37°C in a total reaction volume of 10 μl .

Unit Definition Assay: Two fold dilutions of IdeZ Protease (IgG-specific) are incubated with 1 μg of human IgG and 1X GlycoBuffer 2 in a 10 μl reaction. The reaction mix is incubated for 15 minutes at 37°C . Separation of reaction products are visualized by SDS-PAGE.

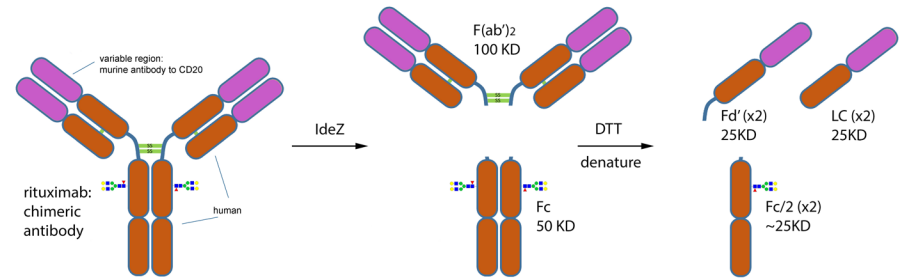

Digestion of IgG with IdeZ Protease (IgG-specific), followed by denaturation.

Molecular Weight: 35,578 daltons

Quality Assurance: No contaminating non-specific proteolytic activity could be detected.

Quality Control Assays

Physical Purity: Purified to > 95% homogeneity as determined by SDS-PAGE analysis using Coomassie Blue detection

Protease Assay: After incubation of 800 units of IdeZ Protease (IgG-specific) with 24 μg of a standard mixture of proteins in a 20 μl reaction, for 20 hours at 37°C , no non-specific proteolytic activity could be detected by SDS-PAGE with Coomassie Blue Detection.

Heat Inactivation: 65°C for 10 minutes.

CERTIFICATE OF ANALYSIS



CERTIFICATE OF ANALYSIS

Reaction Protocol:

The quantity of enzyme recommended is sufficient for the deglycosylation of 50 µg of IgG under non-denaturing conditions. Reactions may be scaled-up linearly to accommodate larger amounts of IgG. Optimal incubation times may vary for a particular substrate.

Typical Reaction Conditions are as Follows:

1. Combine up to 50 µg of human IgG and H₂O in a total reaction volume of 25 µl.
2. Add 2.5 µl of GlycoBuffer 2 (10X).
3. Add 1 µl of IdeZ Protease (IgG-specific).
4. Incubate at 37°C for 30 minutes.*
5. Prepare IgG sample for SDS-PAGE or mass spectrometry analysis.

***Note:** Extend incubation time to 2 hours for murine IgG2a.

Storage: Store at –20°C for up to two years. Avoid multiple freeze/thaw cycles.

Notes:

IdeZ Protease efficiently cleaves human, humanized, chimeric, sheep, rabbit and monkey IgG as well as mouse IgG2a and IgG3. IdeZ Protease will also cleave Fc-fusion proteins, such as Enbrel.

IdeZ Protease does not cleave mouse IgG1 or IgG2b, rat, porcine, bovine or goat IgG. It also does not cleave non-IgG isotypes including IgA, IgM, IgD and IgE.



NEW ENGLAND BIOLABS® is a registered trademark of New England Biolabs, Inc.

This product is intended for research purposes only. This product is not intended to be used for therapeutic or diagnostic purposes in humans or animals.

Companion products

Rapid PNGase F Antibody Standard (P6043)

Rapid PNGase F (P0710)

Remove-iT® Endo S (P0741)

Trypsin-ultra™, Mass Spectrometry Grade (P8101)